





Perform 16/32 samples extraction in approximately 9 mins



AUTOMATED NUCLEIC ACID EXTRACTOR

# MGISP-NE32







Leading Life Science Innovation

MGI Tech Co., Ltd

## ■ MGISP-NE32

MGISP-NE32 is an automated nucleic acid extractor adopting magnetic rod technology. Together with nucleic acid extraction kit, MGISP-NE32 is able to extract and purify nucleic acid from 16/32 samples in approximately 9 mins.

## **■** Technical Principle

The nucleic acid extraction and purification is performed in four automated steps – lysis, binding, washing and elution. Using magnetic rods and disposable tips, MGISP-NE32 is intended for automated moving and processing of magnetic particles in a 96-well plate.



Sample lysis, nucleic acid binding with magnetic particles.



Magnetic particles washing, removing contaminants and inhibitors.



Nucleic acid separation with magnetic particles and elution in elution buffer.

#### 8-inch Built-in Touch Screen

Create, edit and delete program

#### **Reagent and Consumable**

96-well pre-packed plate Disposable tip



 Adjustable mixing module with multiple modes & gears, compatible with different magnetic particles and reagents.

#### **Efficient**

 Moving and processing magnetic particles, extracting and purifying nucleic acid from 16/32 samples in approximately 9 mins.

#### Safe

 UV lamp, 96-well pre-packed plate and disposable tip, avoiding cross-contamination by minimizing contact with sample.

#### **Stable**

 Specially designed heating block for 96-well plate, ensuring even heating during the experiment.



## Temperature Control


### Module

#### **Mixing Module**

Adjustable mixing module with multiple modes & gears



## Performance

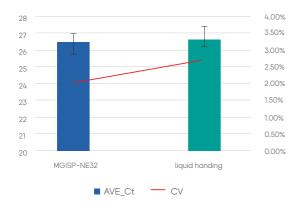

#### Test method

MGISP-NE32 VS Liquid handling method

#### Evaluation standard

RT-PCR

#### Test samples

Simulated samples of throat swab containing novel corona virus (RNA), dilute them in different concentrations

#### Test results

Ct value of fluorescent quantitative PCR is almost equal, that demonstrated the yield of virus RNA is coincident between MGISP-NE32 and liquid handling method. Meanwhile, the reproducibility is great, the CV of Ct value is less than 2.5%.

## ONE-STOP SOLUTION

Provide instrument, consumables and kits for extraction



#### **MGISP-NE32**

- Automated Nucleic Acid Extractor
- Automated extraction step
- Moving and processing magnetic particles
- Nucleic acid extraction from 16/32 samples in approximately 9 mins.



#### Nucleic Acid Extraction Kit

- 96-well pre-packed plate
- Good extraction stability and reproducibility
- Compatible with Automated Nucleic Acid Extractor
- Adapted to most downstream applications:
  PCR, RT-PCR, Sequencing, etc.

# **MGISP-NE32**

## Automated Nucleic Acid Extractor

Perform 16/32 samples extraction in approximately 9 mins

#### **ORDER INFORMATION**

#### **Equipment**

Cat. No.	Product name	Specification
950-000020-00	MGISP-NE32RS Automated Nucleic Acid Extractor	EA, CE RUO
950-000013-00	MGISP-NE32 Automated Nucleic Acid Extractor	EA, CE IVD

#### Reagent

Cat. No.	Product name	Specification	Feature
1000023774	MGIEasy Nucleic Acid Extraction Kit (OP02-32)	32 preps, CE RUO	35 mins, Standard Version
1000022606	Nucleic Acid Extraction Kit (OP02-32)	32 preps, CE IVD	35 mins, Standard Version
1000023938	MGIEasy Nucleic Acid Extraction Kit (VDR03P-32)	32 preps, CE RUO	9 mins, Fast Version
1000023937	Nucleic Acid Extraction Kit (VDR03P-32)	32 preps, CE IVD	9 mins, Fast Version

#### Copyright Disclaimer

The copyright of this brochure is solely owned by MGI Tech Co. Ltd.. The information included in this brochure or part of, including but not limited to interior design, cover design and icons, is strictly forbidden to be reproduced or transmitted in any form, by any means (e.g. electronic, photocopying, recording, translating or otherwise) without the prior written permission by MGI Tech Co., Ltd.. All the trademarks or icons in the brochure are the intellectual property of MGI Tech Co., Ltd. and their respective producers.

Version: December 2020 | MGPC4002202



https://bit.ly/34Cfqpe